




## **Removal of Compressor Shelf**

From  
Computerized Environmental Chamber

**NOTE: DISCONNECT CHAMBER FROM POWER SOURCE**

1. Remove top from chamber.
2. Remove right-side top panel by removing the screw in the bottom flange (Grounding Screw) then lifting upwards and pulling outwards.
3. Disconnect Power Lead to Compressor Shelf. (3 Wires located at the Terminal Block, Color Coded: Green-White-Black.) See Drawing.
4. Disconnect Heated Control Wire. (At Quick Disconnect.)
5. Remove the three (3) screws located on the bottom side of Compressor Shelf.
6. The Complete Compressor Shelf can now be slid out.

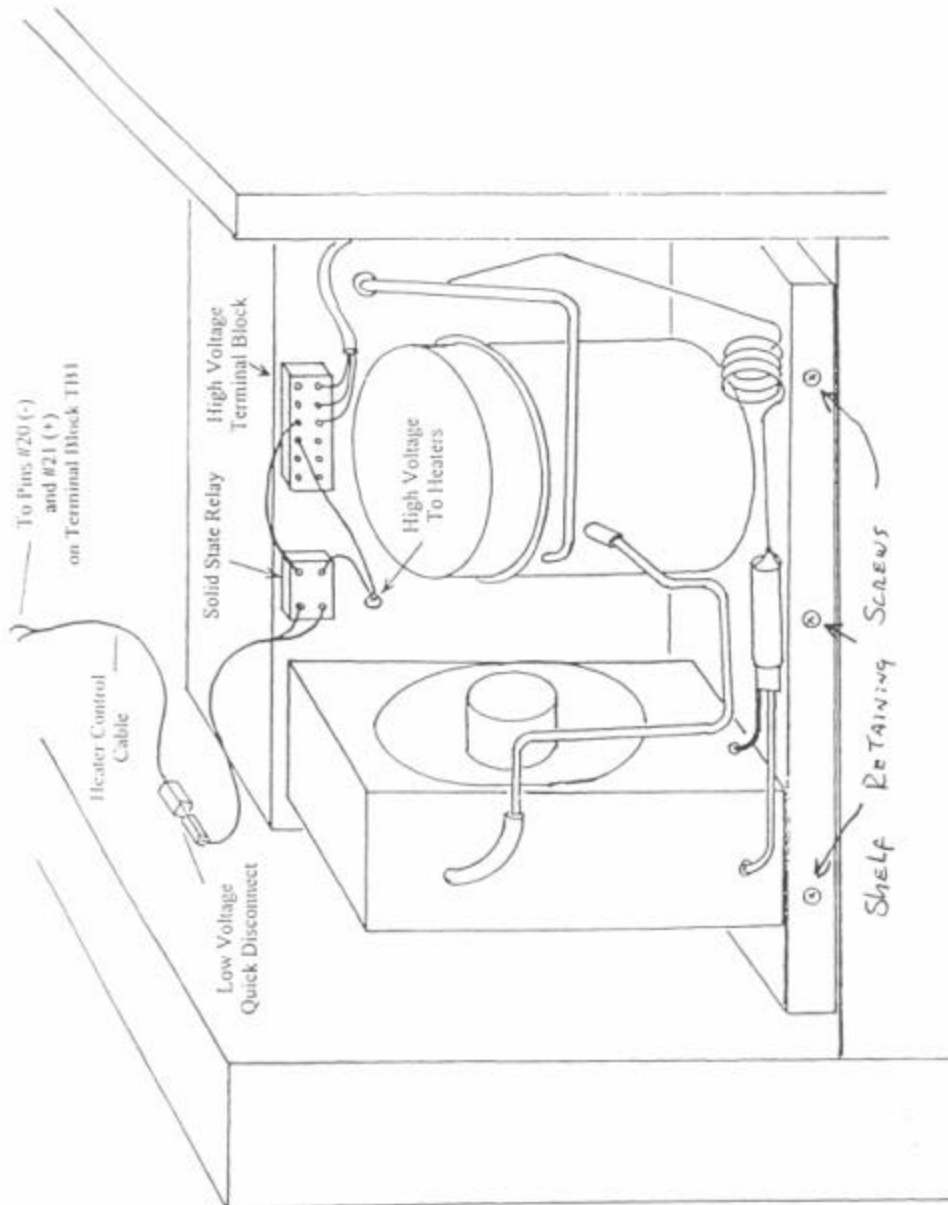

**Innovative Design, Engineering and Manufacturing since 1946**

4420 San Fernando Road Glendale CA 91204 818 247 7500 fax 818 247 7257 melco@melcowire.com www.melcowire.com